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Gamma - SIM

File MCA Calibrate Display Analyze Edit Options Datasource Help

Acquire Start F4  
Acquire Stop F4  
Acquire Setup...  
Clear  
**Adjust...**  
Status  
Advance  
Load...

1.9 keV Counts: 0 Preset: 3600/134.31 VFS-64

Energy(keV)

TIME INFO

	Acq. Start:	2003-4-21 14:09:51	Elapsed	Preset
Next	Dead Time:	0.51%	Live (secs.): 134.31	3600
Prev	Comp. Preset Region:	0 - 0 (channels)	Real (secs.): 135.00	0
			Total (cnts.): 0.00	0

Adjust device settings Execution Status: ready

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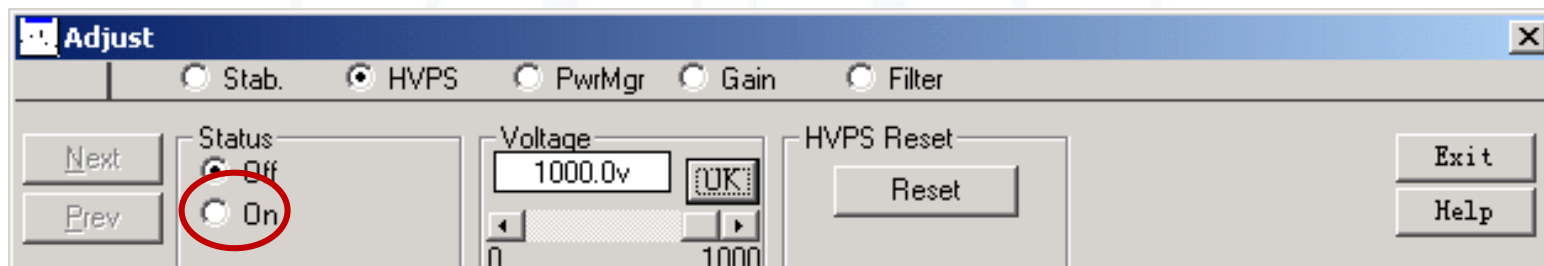
选择Analyze中Peak Locate的Unidentified 2nd Diff...

TIME INFO

	Acq. Start:	Live (secs.):	Elapsed	Preset
Next	Dead Time: 0.52%	44.768	44.768	1000
Prev	Comp. Preset Region:	Real (secs.): 45.000	0.00	0
	0 - 0 (channels)	Total (cnts.):		0

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Peak Locate Unidentified 2nd Diff. Setup

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选择Analyze中Peak Area的Sum/Non-Linear LSQ Fit...

TIME INFO

	Acq. Start:	Elapsed	Preset
Next	Dead Time: 0.52%	Live (secs.): 44.768	1000
Prev	Comp. Preset Region:	Real (secs.): 45.000	0
	0 - 0 (channels)	Total (cnts.): 0.00	0

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**Sum / Non-Linear LSQ Fit Setup**

Peak Area Region  
Start channel:   
Stop channel:

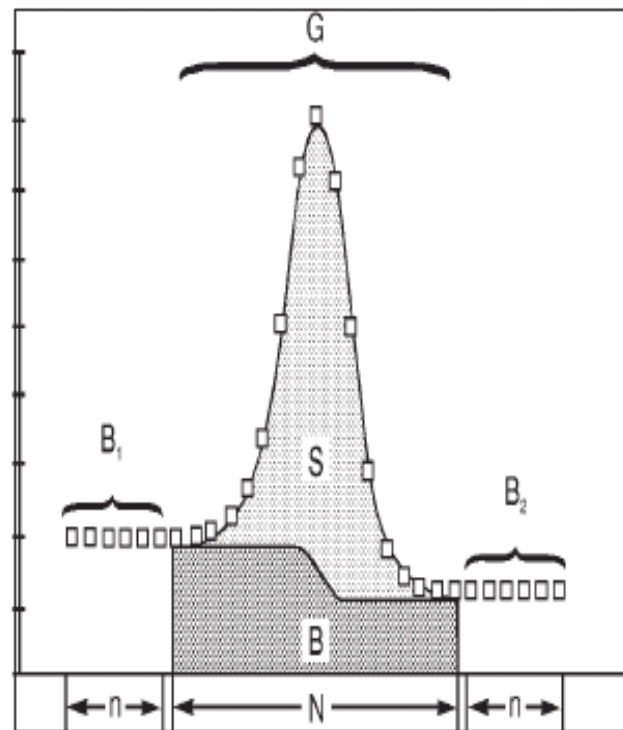
Continuum:  Channels  
 Channels  FWHM

Continuum function:

Residual Search  
 Perform Search    Threshold:   
Minimum separation (FWHM):

ROI Limits Determination  
Max. Num. FWHMs between peaks:   
Max. Num. FWHMs for left limit:   
Max. Num. FWHMs for right limit:

95% Critical level test  
 Use fixed FWHM  
 Use fixed tail parameter  
 Fit singlets  
 Display ROIs  
 Reject zero area peaks

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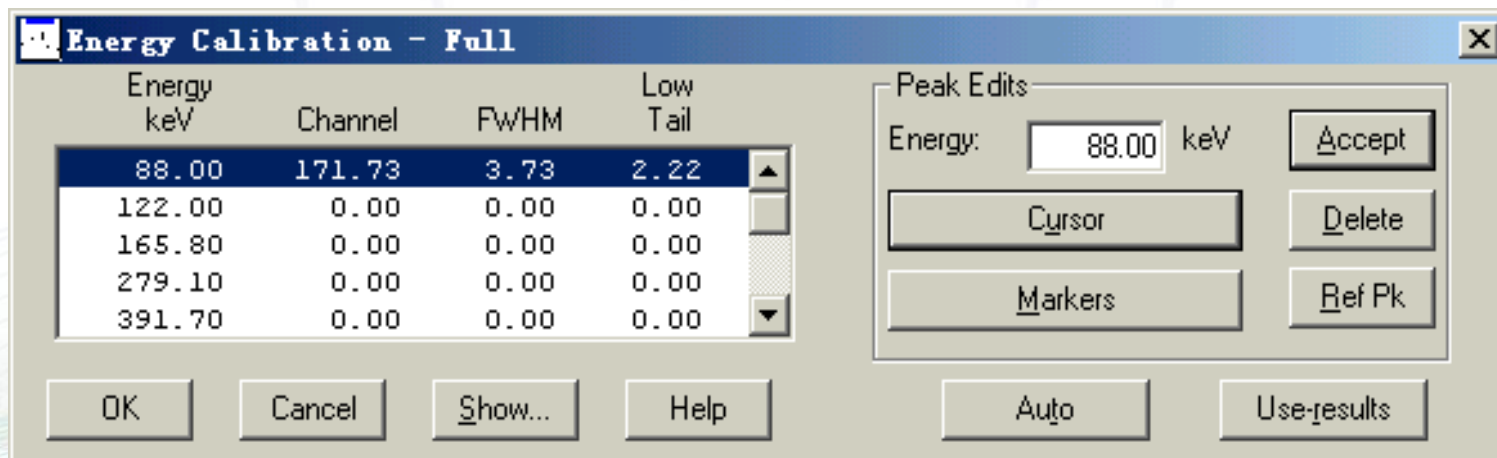
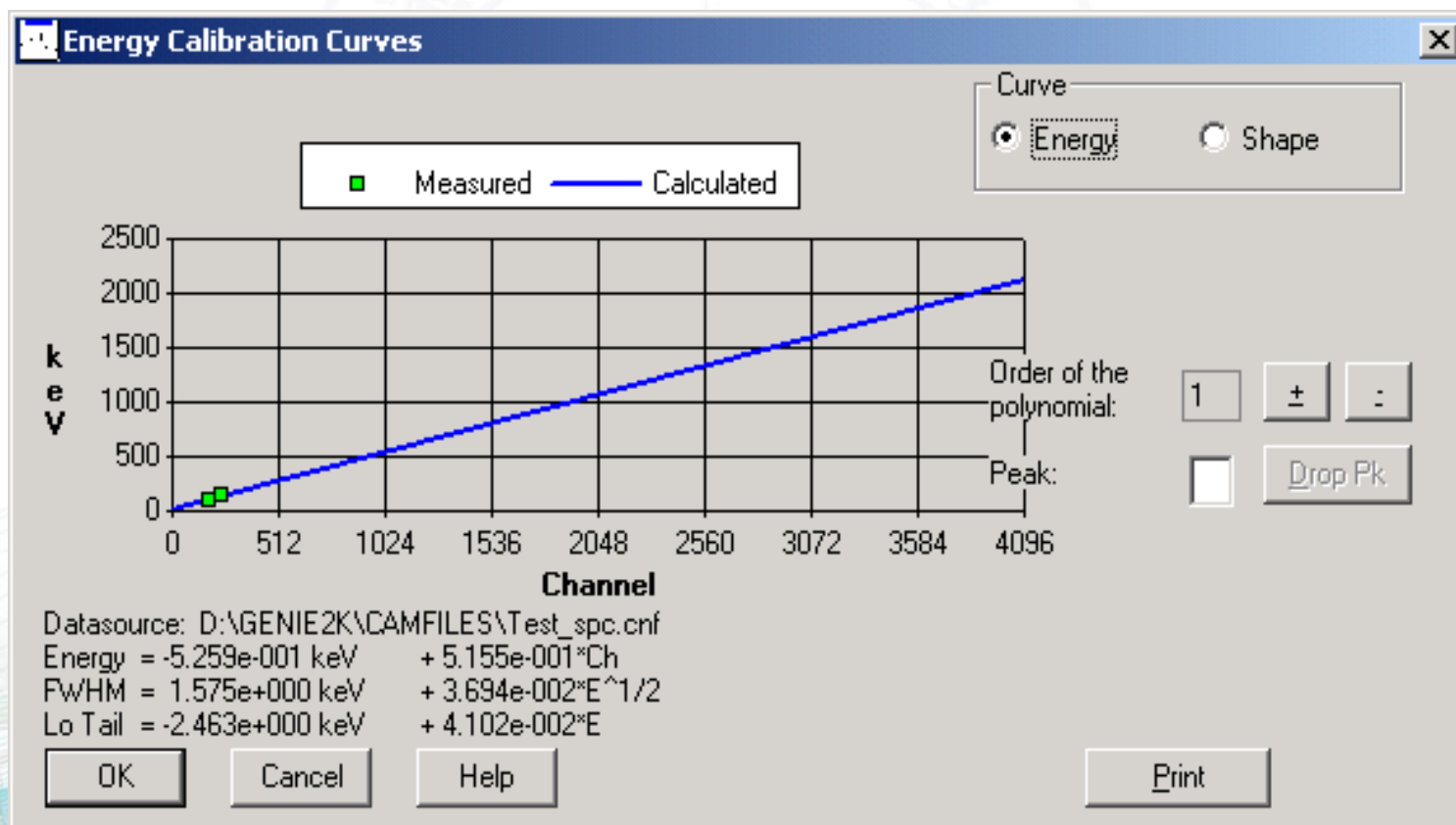
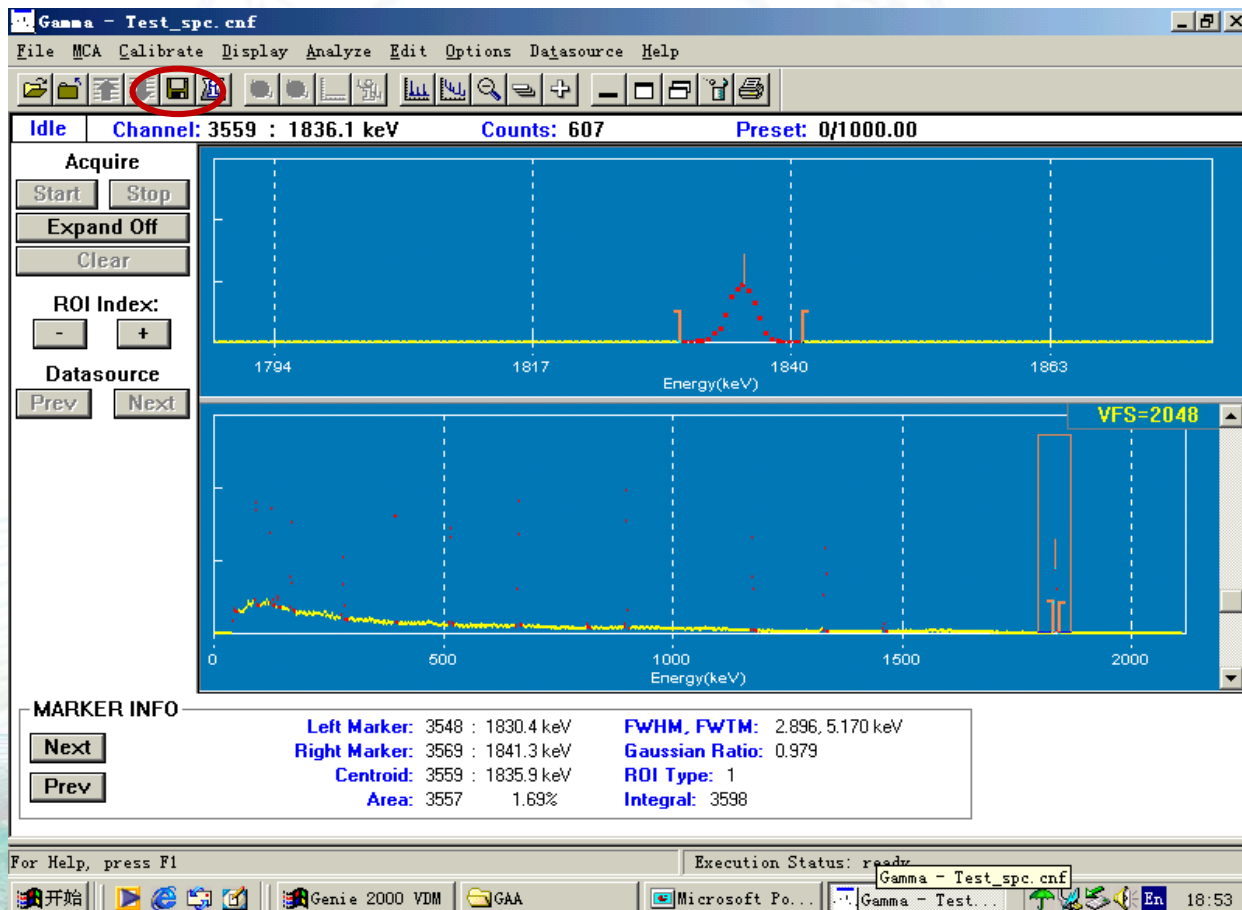
选择Calibrate中Efficiency的By Certificate File

The screenshot shows the Gamma software interface. The 'Calibrate' menu is open, and 'Efficiency' is selected. The 'By Certificate File...' option is highlighted with a red circle. The main window displays a gamma spectrum with markers at 1840 and 1863 keV. The 'MARKER INFO' table is visible at the bottom.

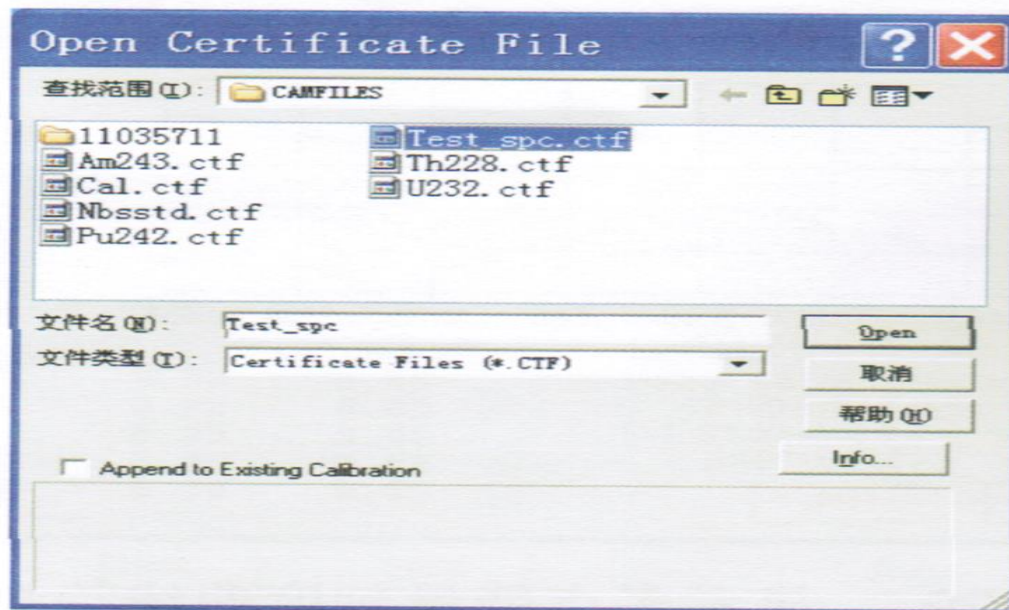
MARKER INFO	
Left Marker:	3548 : 1830.4 keV
Right Marker:	3569 : 1841.3 keV
Centroid:	3559 : 1835.9 keV
Area:	3557 : 1.69%
FWHM, FWTM:	2.896, 5.170 keV
Gaussian Ratio:	0.979
ROI Type:	1
Integral:	3598

Execution Status: ready

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启动Calibrate/Efficiency/By Certificate File菜单命令，出现以下界面

Calibrate/Efficiency/By Certificate File对话框

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
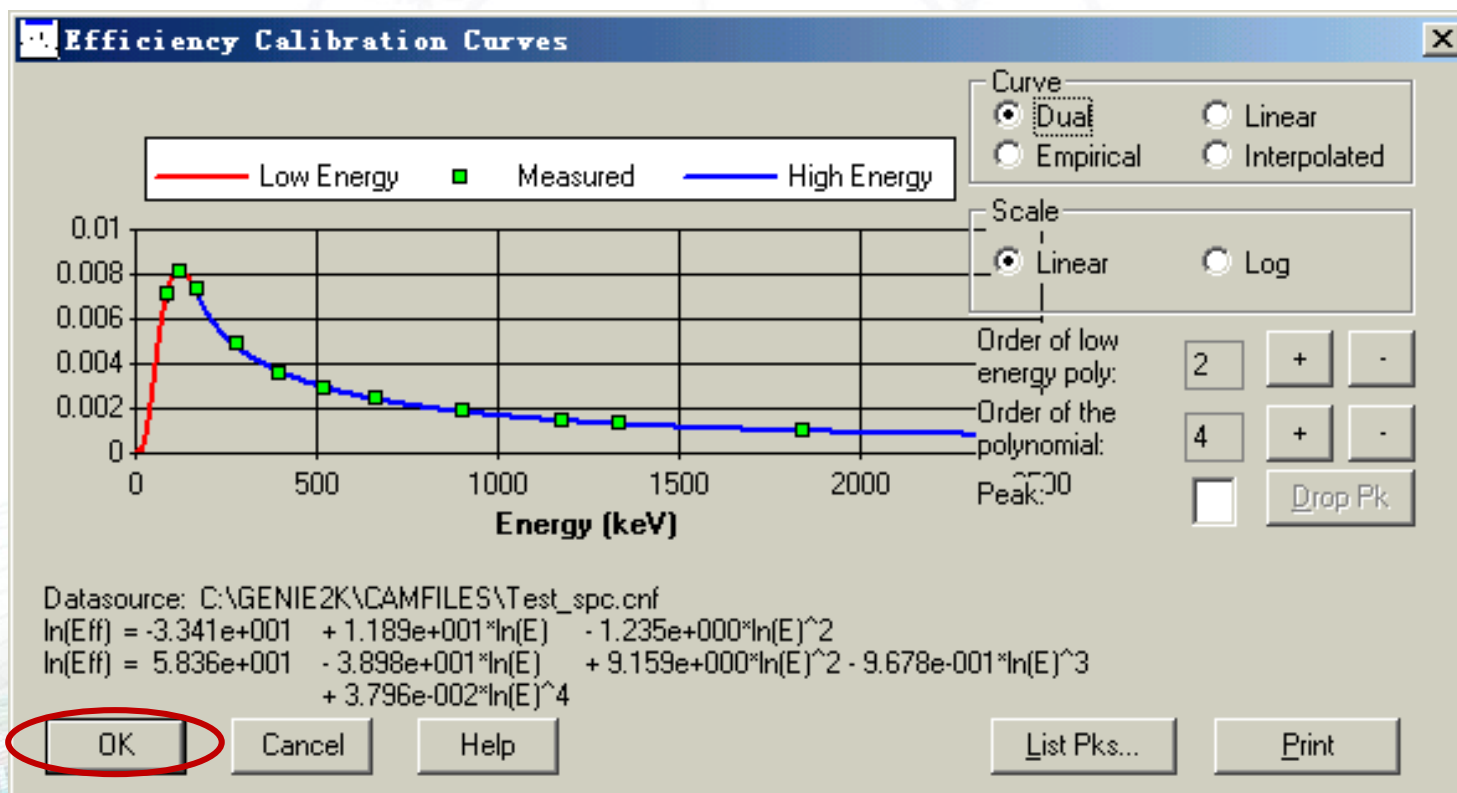
The screenshot shows the Gamma software interface. The 'Calibrate' menu is open, and the 'Store...' option is highlighted with a red circle. The main window displays a gamma spectrum plot with peaks and a 'TIME INFO' table at the bottom.

TIME INFO	Acq. Start:	1983-5-5 8:53:00	Elapsed	Preset
Next	Dead Time:	0.50%	Live (secs.):	1000.000
Prev	Comp. Preset Region:	0 - 0 (channels)	Real (secs.):	1005.000
			Total (cnts.):	0.00

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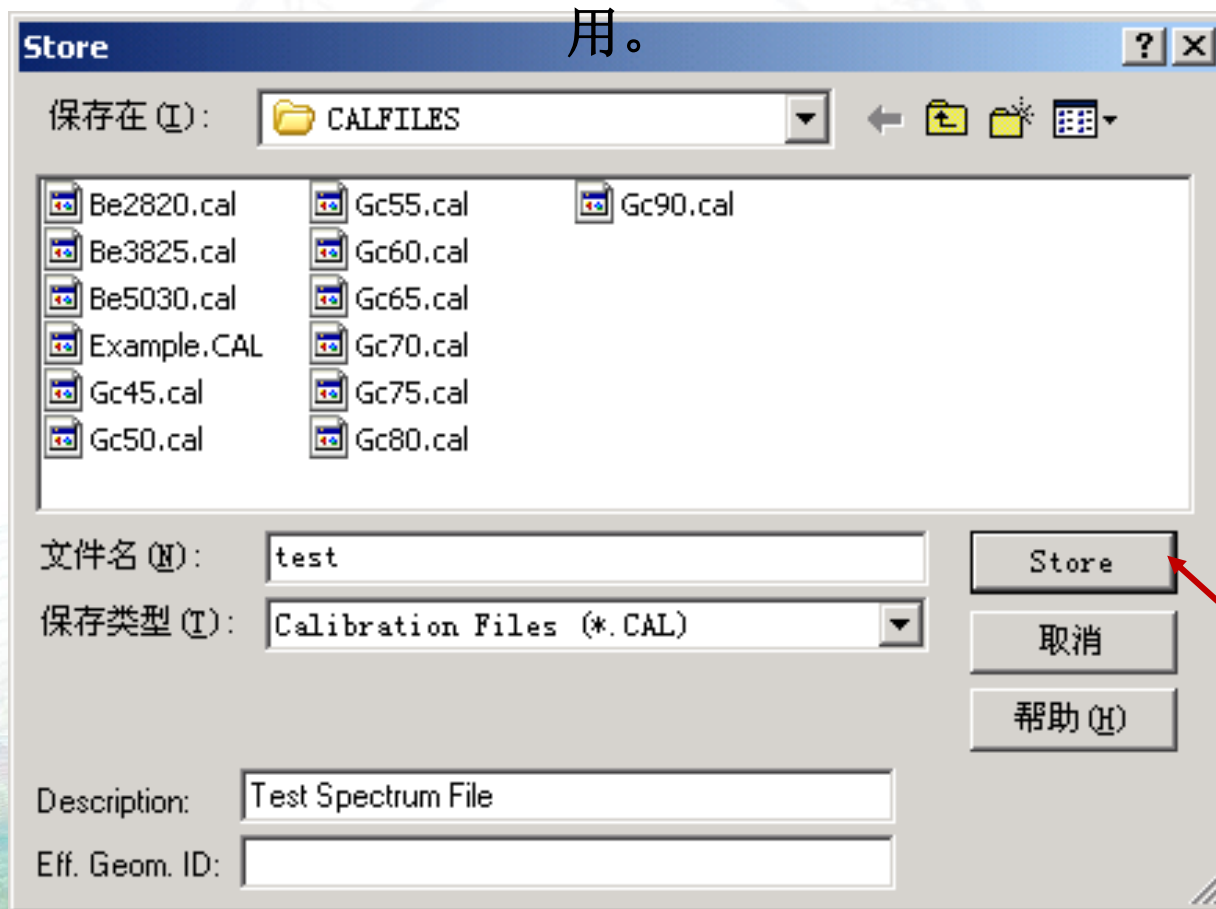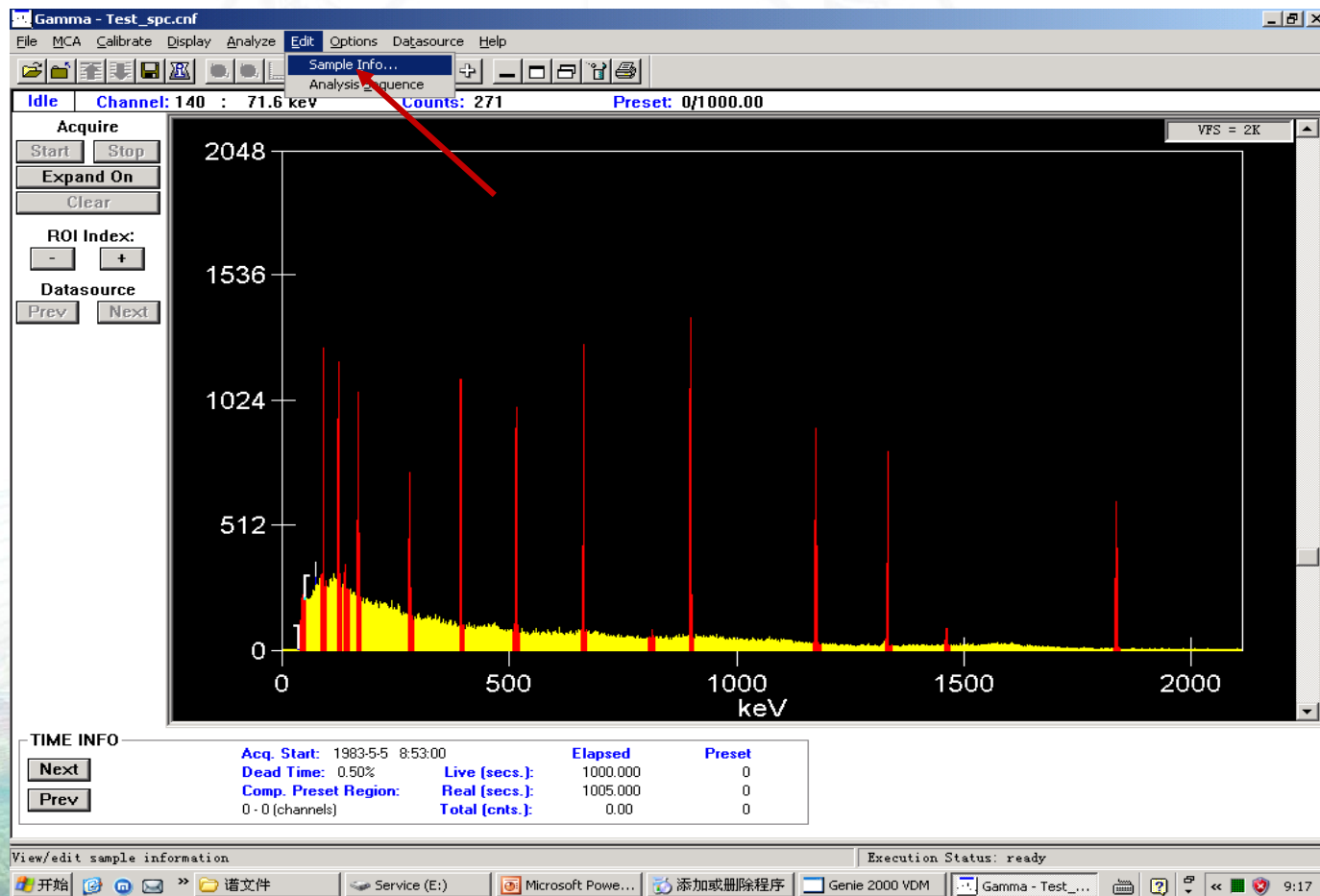
**Edit Sample Information**

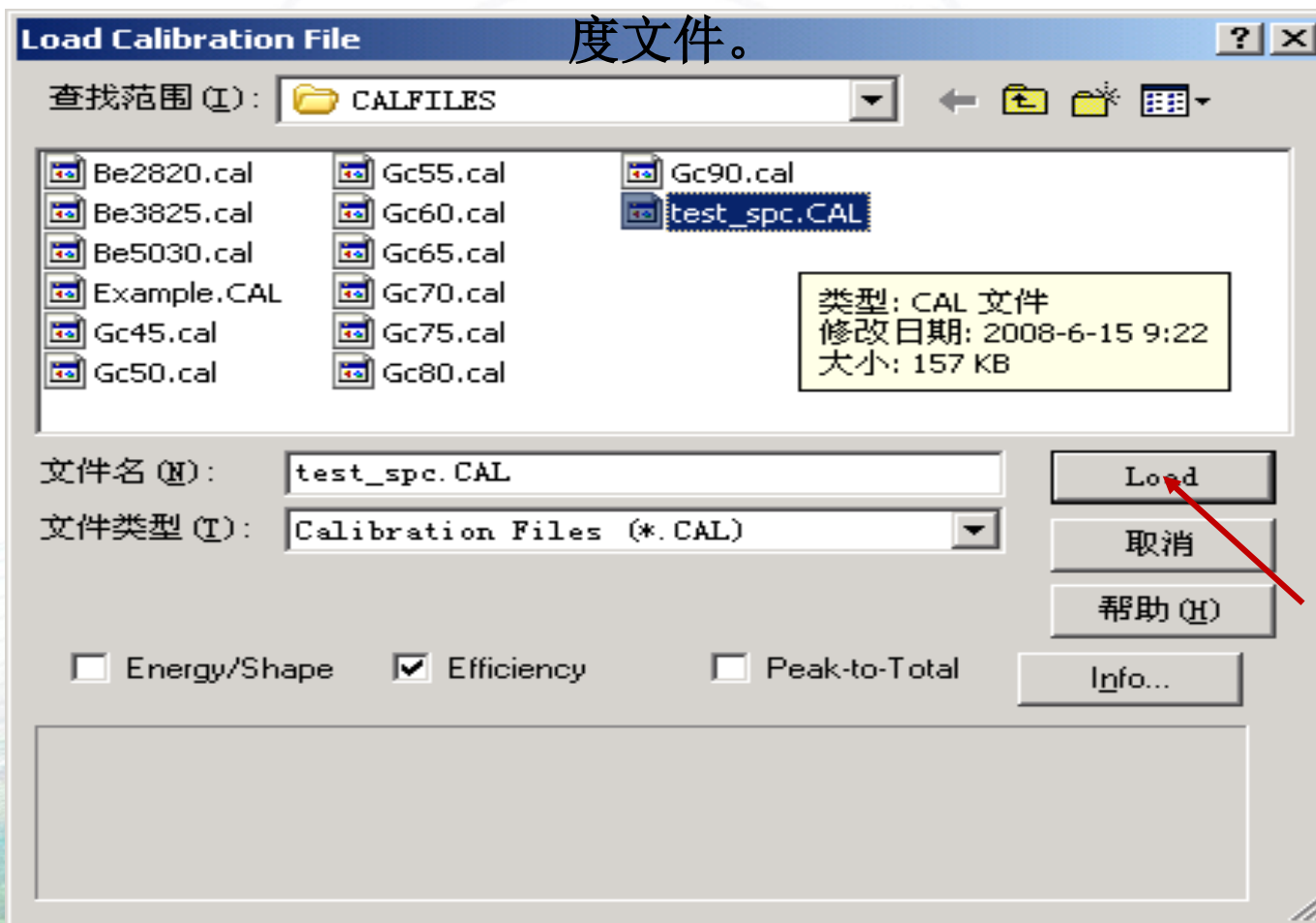
Sample Title:	<input type="text" value="GENIE-PC Spectrum No. 2"/>	Sample ID:	<input type="text" value="0629QB #1"/>
Collector Name:	<input type="text" value="M. J. Koskelo"/>	Type:	<input type="text"/>
Sample Description:	<input type="text" value="Derived from APOGEE File 505"/> <input type="text"/> <input type="text"/>	Quantity:	<input type="text" value="1"/>
Buildup Type:	<input checked="" type="radio"/> None <input type="radio"/> Deposition <input type="radio"/> Irradiation	Uncertainty:	<input type="text" value="0"/>
Begin Date:	<input type="text"/> at <input type="text"/>	Units:	<input type="text" value="Unit"/>
Sample Date:	<input type="text" value="1983-2-1"/> at <input type="text" value="8:00:00"/>	Sample Geometry:	<input type="text" value="Geometry #1"/>
		Random Error (%):	<input type="text" value="0"/>
		Systematic Error (%):	<input type="text" value="0"/>

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选择File中的Save As

The screenshot shows the Gamma software interface. The 'File' menu is open, and 'Save As...' is highlighted with a red circle. The main window displays a gamma spectrum plot with a peak at 1840 keV. The plot shows Energy (keV) on the x-axis and Counts on the y-axis. A Gaussian fit is overlaid on the peak. The 'MARKER INFO' section at the bottom provides detailed data for the peak.

MARKER INFO	
Left Marker:	3548 : 1830.4 keV
Right Marker:	3569 : 1841.3 keV
Centroid:	3559 : 1835.9 keV
Area:	3557 : 1.69%
FWHM, FWTM:	2.896, 5.170 keV
Gaussian Ratio:	0.979
ROI Type:	1
Integral:	3598

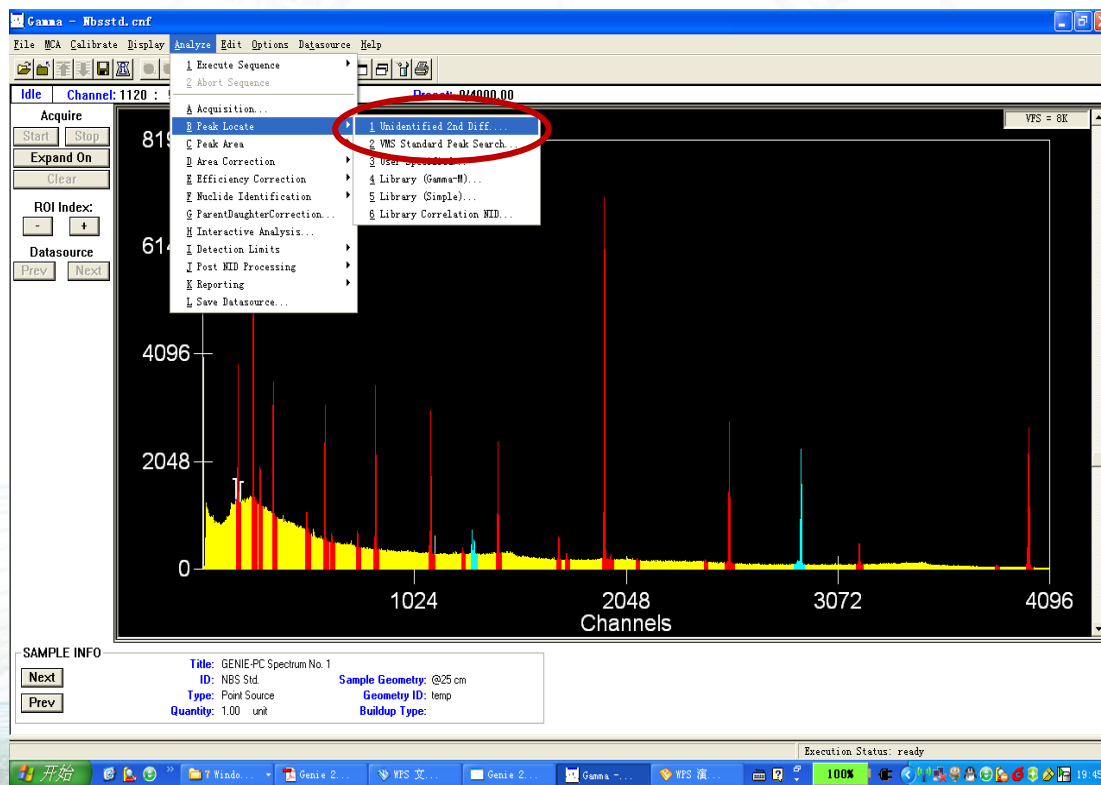
Execution Status: ready

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寻峰(Peak Locate)推荐参数—Unidentified 2nd Diff

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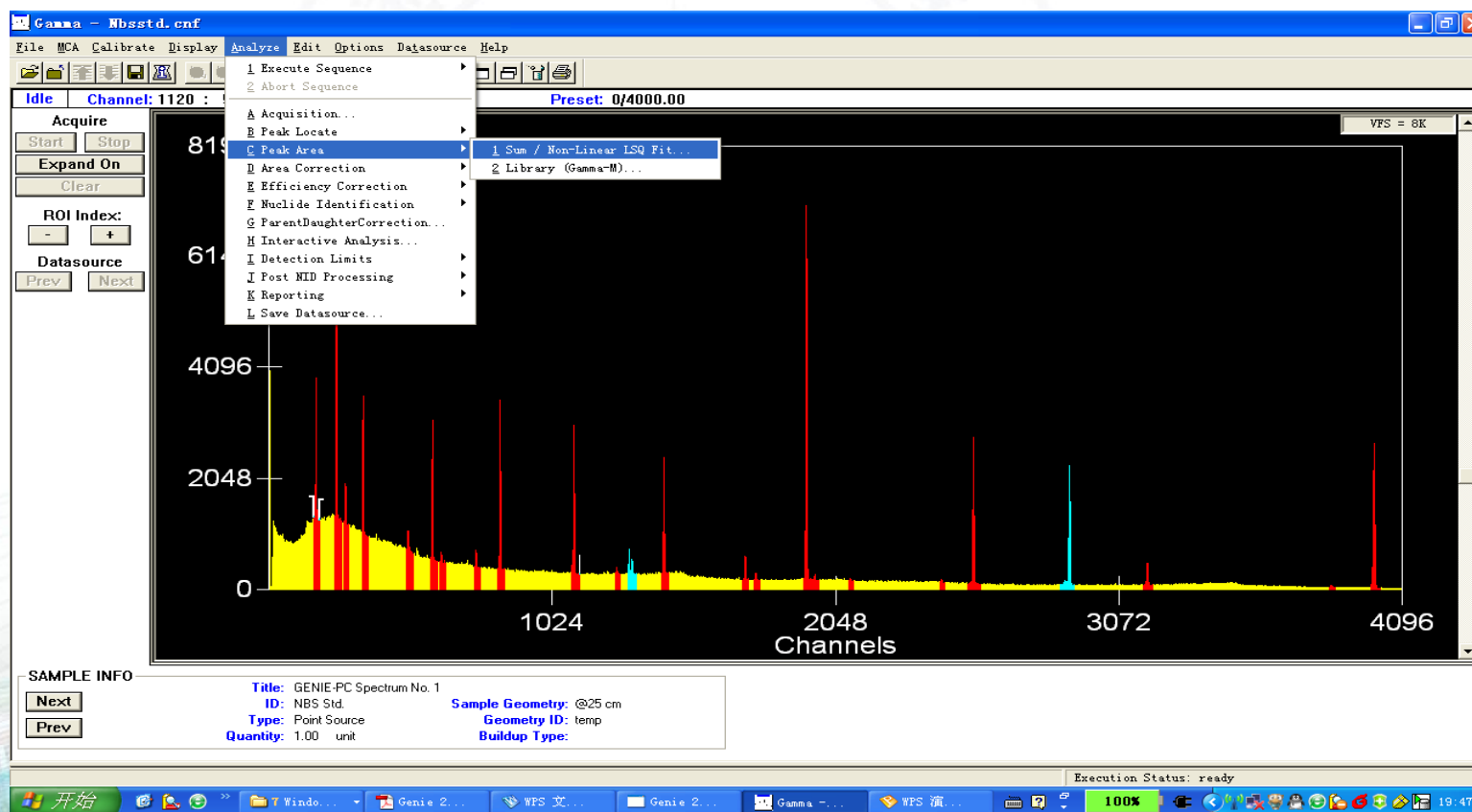
峰面积(Peak Area)推荐参数—Sum/Non-Linear Least Squares Fit Peak Area

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峰面积(Peak Area)推荐参数—Sum/Non-Linear Least Squares Fit Peak Area

**Sum / Non-Linear LSQ Fit Setup**

Peak Area Region  
Start channel:   
Stop channel:

Continuum:  Channels  
 Channels  FWHM

Continuum function:

Residual Search  
 Perform Search Threshold:   
Minimum separation (FWHM):

ROI Limits Determination  
Max. Num. FWHMs between peaks:   
Max. Num. FWHMs for left limit:   
Max. Num. FWHMs for right limit:

95% Critical level test  
 Use fixed FWHM  
 Use fixed tail parameter  
 Fit singlets  
 Display ROIs  
 Reject zero area peaks

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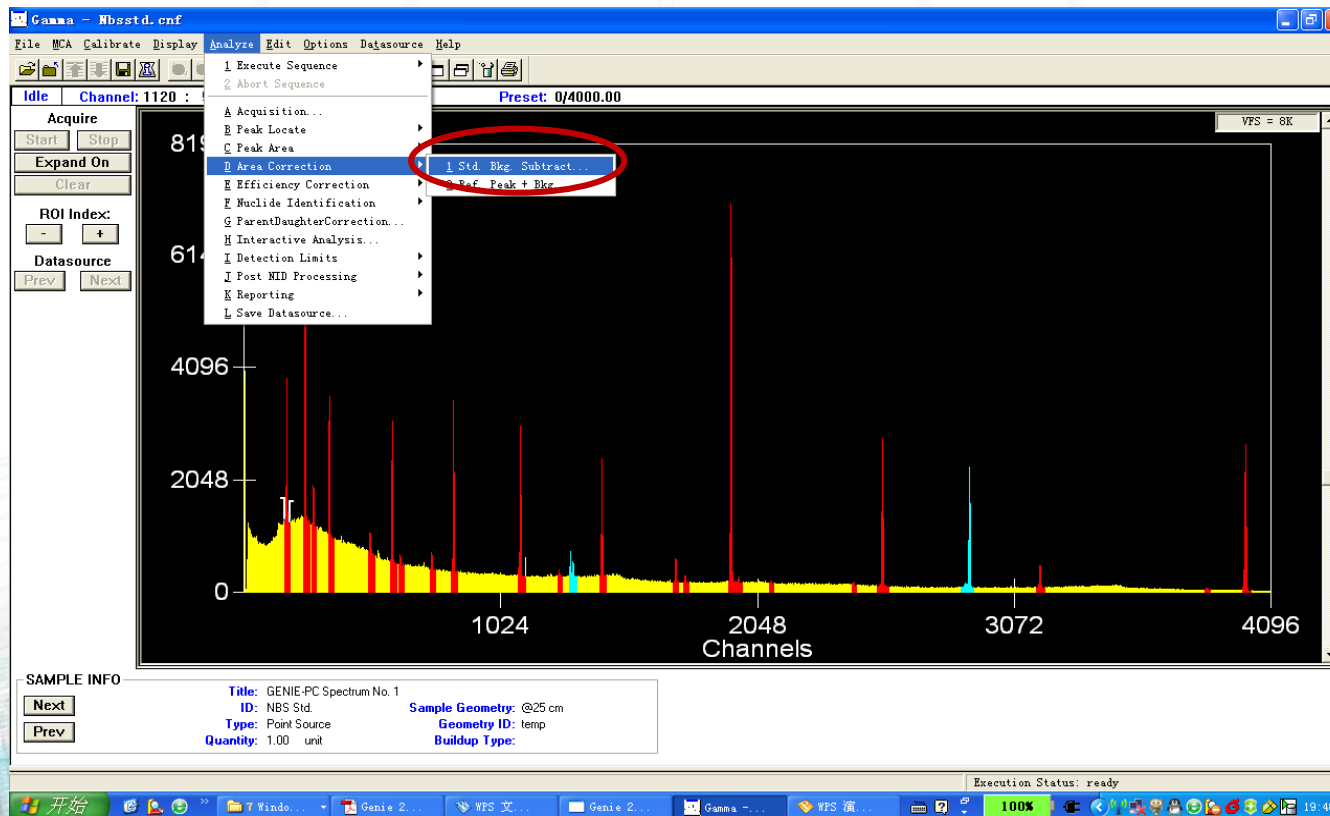
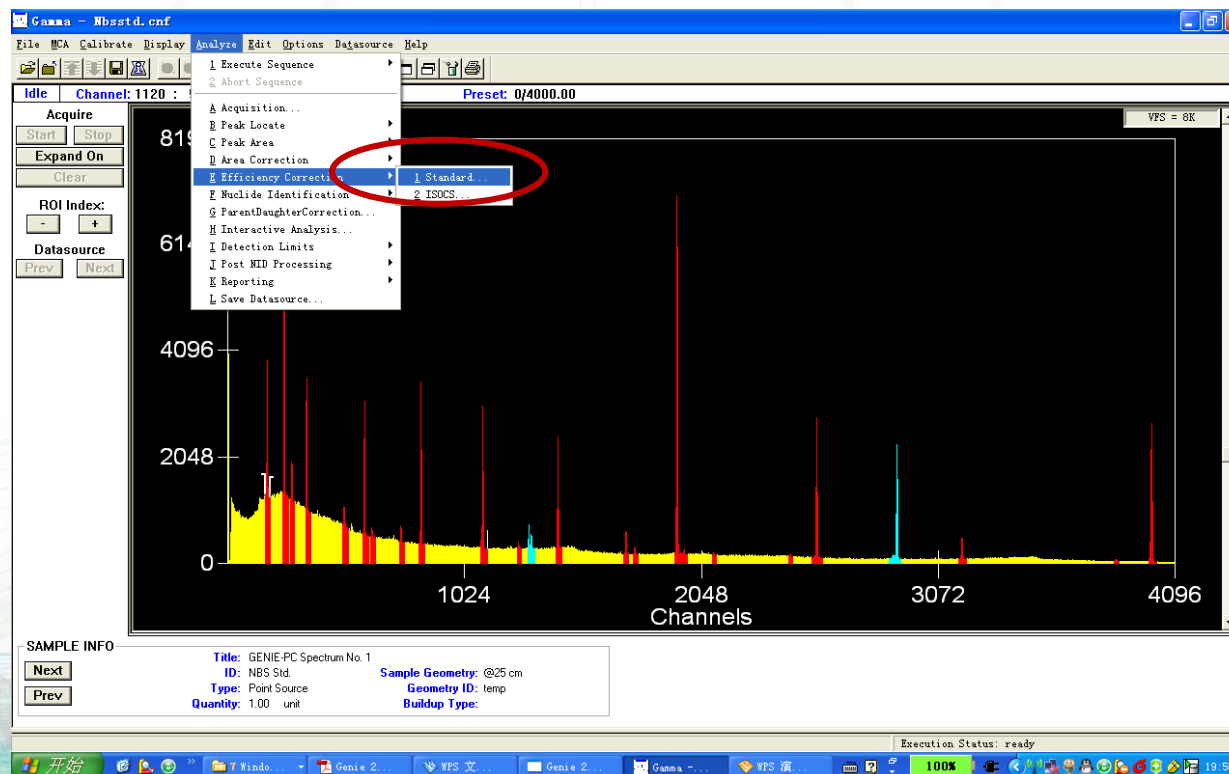
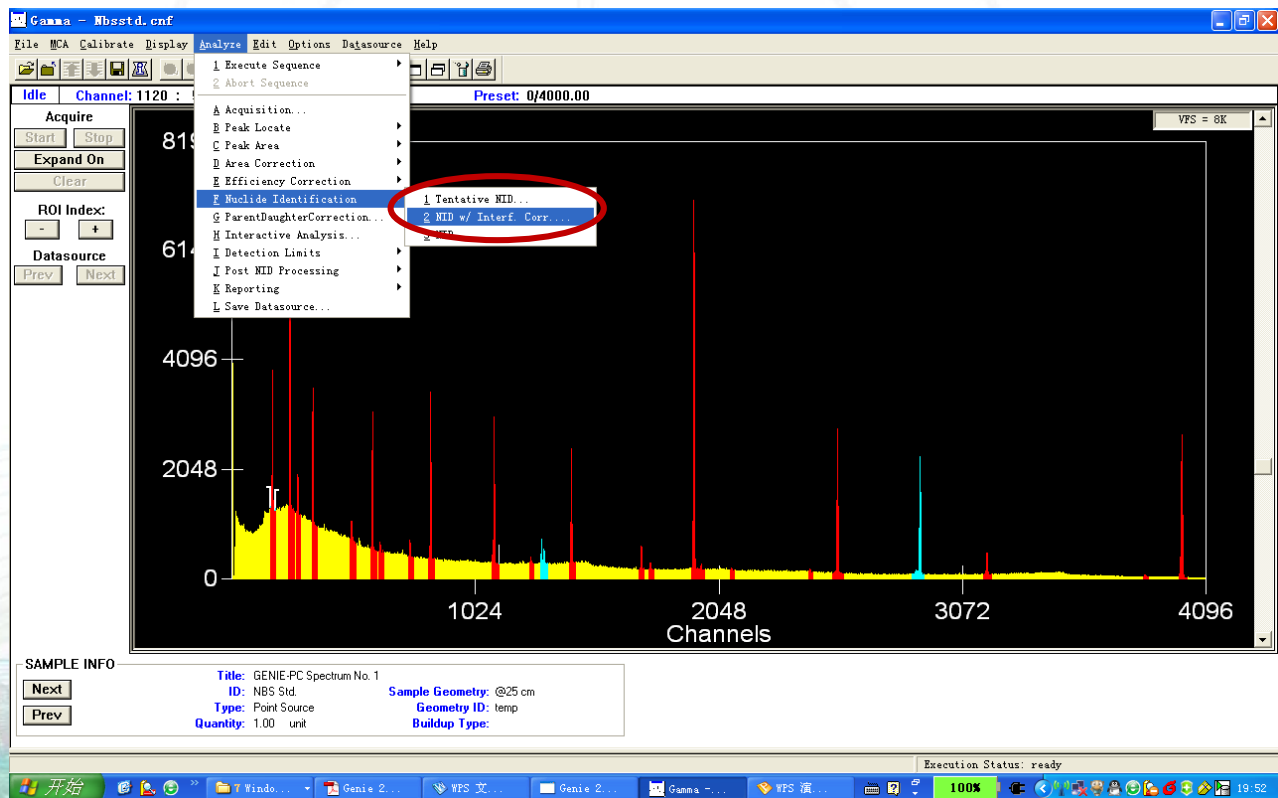
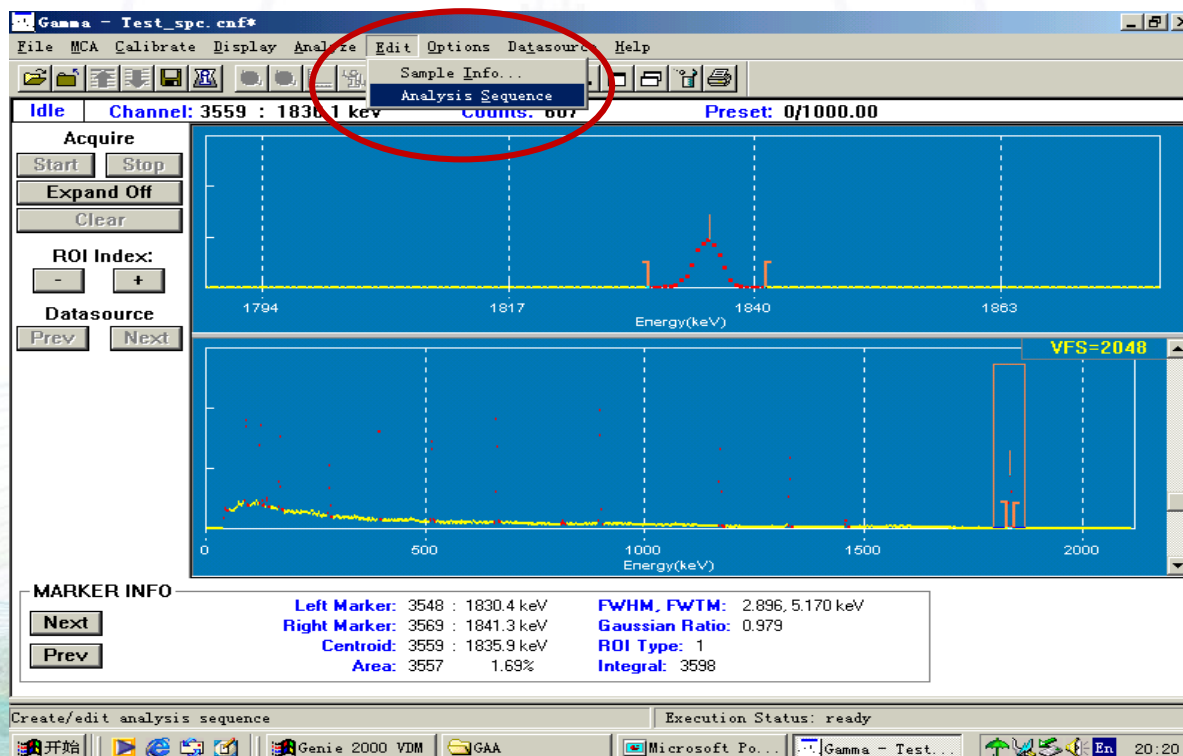
从Analyze的Execute Sequence中可随时调用

## Basic Analysis

The screenshot shows the Gamma software interface. The 'Analyze' menu is open, and 'Basic Analysis' is highlighted. The main window displays a gamma spectrum with a peak at 40.7 keV. The 'MARKER INFO' section shows the following data:

Parameter	Value
Left Marker	71 : 35.9 keV
Right Marker	84 : 42.6 keV
Centroid	80 : 40.7 keV
Area	312 : 15.42%
FWHM, FWTM	0.661, 4.229 keV
Gaussian Ratio	3.510
ROI Type	1
Integral	1705

The 'MARKER INFO' section also includes 'Next' and 'Prev' buttons. The 'Execution Status' is 'ready'. The taskbar shows the system time as 23:45.